Kit contains:
4 sterile test vials
Test Instructions

24-Hour In-Office Dental Waterline Test Kit

Detects microbial levels that exceed 500 CFU/mL
Easy-to-Read Color Change Results

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REF **IMS-1004**

(01)10889950140014

HuFriedyGroup
The Best In Practice

If you are interested in learning more about the
GreenLight Dental Compliance Center™ by
Hu-Friedy to track your in-office dental waterline test
results, scan the QR code.

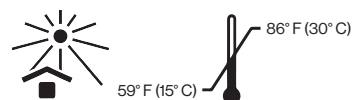
Hu-Friedy

Manufactured for:
Hu-Friedy Mfg. Co., LLC
2501 Sanders Rd.
Northbrook, IL 60062 | USA
1-800-Hu-Friedy

Made in Canada

N5568 / REV 4 - 2512

All company and product names are trademarks of Hu-Friedy Mfg. Co., LLC, its affiliates or related companies, unless otherwise noted.
©2025 Hu-Friedy Mfg. Co., LLC. All rights reserved.

Directions for Use

Scan the QR code for support with the aseptic technique for obtaining a
water sample for testing.

- Using an indelible ink marker, number the cap(s) of the vial(s) selected for testing and log* that number along with the name of the clinician drawing the sample, sampling location (i.e.: Op 1, Hygiene Room etc.), source (i.e.: AWS, high speed handpiece etc.), date and time water sample was taken. For Example:

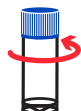
- Vial 1: M. Summers, Op 1, AWS, 3/15/23, 9:30am
- Vial 2: M. Summers, Op 2, AWS, Scaler, High Speed Handpiece, 3/15/23, 9:40am

* Scan the QR code to download a complimentary Waterline Testing log.

- Perform hand hygiene and don disposable treatment gloves.

- Remove the cap from the first vial to be used for sample collection. Keep the cap and vial in your hand, or place the cap open side up on a clean surface.

• To avoid contamination, do not open more than one vial at a time.

Fill selected vials to the 5mL fill line with water from the line(s) to be tested and seal immediately. Do not underfill. *If pooling multiple waterlines into a single vial*, note that the pooled sample:

- Must use equal amounts of water from each waterline
- Include no more than 10 lines
- Pull from a single unit, in a single operatory

4.

Once the vial(s) are sealed, shake well to dissolve the dry media deposited at the bottom of the vial. The water inside the vial will turn blue.

Note: If the water does not become blue after shaking, void the test and contact HuFriedyGroup Customer Care.

It is recommended to take a picture of your test now in front of a sheet of white paper, and note the date and time.

- Incubate the vial(s) at room temperature (no incubator needed) in an upright position for 24-hours.

- Read results no later than 24-hours following sampling.**

- Check for water color change and compare against the color chart below.

● Clean water
0 - 500 CFU/mL

Tip: Holding the test vial up against a sheet of white paper can help in determining color changes.

- Any change from the initial blue water color indicates a microbial level in the water sample above 500 CFU/mL.**

- If a color separation occurs after 24-hours, shake to determine final color results.
- Waterlines with test results above 500 CFU/mL require remediation. For assistance, scan the QR code for HuFriedyGroup failed waterline testing remediation recommendations, contact HuFriedyGroup Customer Care, or email watersupport@hu-friedy.com

It is recommended to take a picture of your results in front of a sheet of white paper at the 24-hour mark (in the same location where you took the initial picture), and attach them to your waterline testing log along with the picture taken when incubation started.

- File and retain these test results for the length of time required by your state dental board.

Test Disposal:

Test with passing results:

- The contents in test vials with passing results may be flushed down any sink drain. The cap and vial may be presented for recycling or discarded in your regular office trash.

Failed Tests:

- It is recommended to autoclave failed tests. Following autoclaving, contents may be flushed down any sink drain. The cap and vial may be presented for recycling or discarded in your regular office trash.

For failed test remediation support, or questions related to dental waterline testing or treatment, please contact Customer Care (800) 483-7433, or email watersupport@Hu-Friedy.com

HuFriedyGroup
The Best In Practice

24-Hour In-Office Dental Waterline Test Kit

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Detects microbial levels that exceed
500 CFU/mL
Easy-to-Read Color Change Results

Kit contains:
4 sterile test vials
Test Instructions

HuFriedyGroup
The Best In Practice

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